



Using IDEXX VetLab* with Cornerstone* Software


This guide explains the recommended way to create an IDEXX VetLab* in-clinic laboratory request and view the test results.

IMPORTANT: This document applies to Cornerstone* 8.4 NEXT or later. Some features may have changed from previous versions. Refer to additional online resources and Cornerstone help  for more information.

Create a laboratory request using the Patient Visit List or an invoice

Note: The invoice item must be linked to the profile, and the **Use special actions** check box must be selected in the INCLINIC laboratory preferences. For details, see the *IDEXX VetLab* with Cornerstone* Software Setup Guide*.

1. On the **Patient Clipboard*** , in the **Patient List** area, right-click the patient name, and then select either **Patient Visit List (PVL)** or **Invoice**.
2. Enter the invoice item ID for the test, and then press **TAB** to open a new laboratory request.

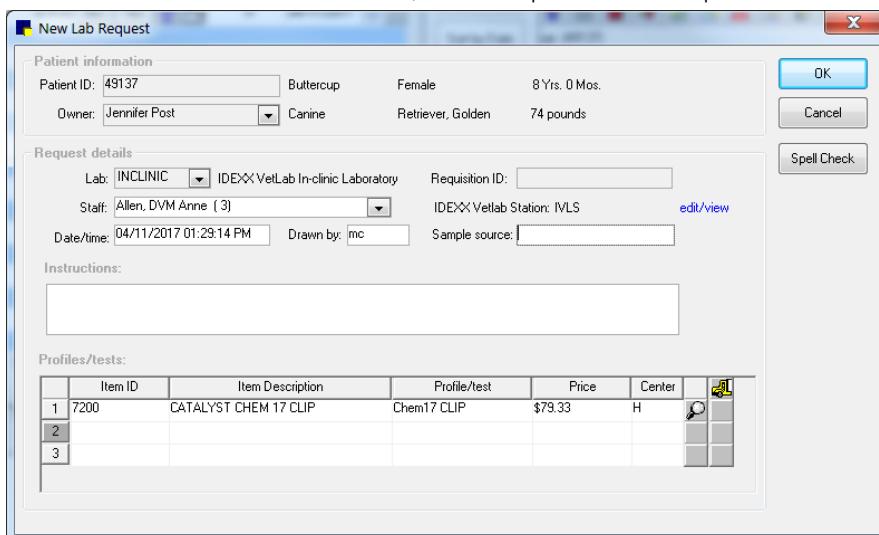

Item ID	Item Description	Profile/test	Price	Center
1 7200	CATALYST CHEM 17 CLIP	Chem17 CLIP	\$79.33	H
2				
3				

3. Enter or modify the staff information as needed.
4. Enter values in the **Drawn by** box, if required, and then click **OK**.
5. If more tests are needed, click **Close**, and then add the additional invoice items.
6. After all tests are visible in the last request, click **OK**.


Running tests on the IDEXX VetLab Station

For detailed information about using the IDEXX VetLab Station, see the *IDEXX VetLab* Station Operator's Guide*.

1. On the IDEXX VetLab Station Home screen, select the patient on the **Pending** tab.
2. Tap **Run**.
3. Confirm the reason for testing.
4. Tap the appropriate instrument icon(s).
5. Tap **Run**.

View results in the patient record using VetConnect* PLUS

If VetConnect* PLUS has been activated, the VetConnect PLUS tab will be available whenever you view laboratory results.

1. On the toolbar, click **Patient Clipboard** .
2. Select the client and patient record.
3. Select the **VetConnect PLUS** tab.

Note: Internet Explorer* 9 or higher is required to view VetConnect PLUS results.


Additional resources

Refer to the following for more information:

- IDEXX Learning Center
- Web Links > Cornerstone
- Cornerstone F1 Help

Find this guide

Look under “Resources by Topic” on the [Cornerstone Software Resources](#) page.

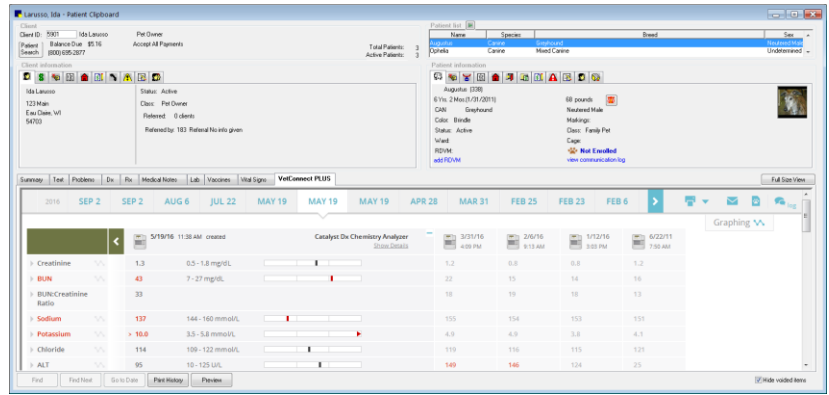

The screenshot displays the IDEXX VetConnect PLUS software interface. At the top, there is a 'Patient Clipboard' window showing patient information for 'Ida Linares', including her ID, balance due, and contact details. Below this, there are tabs for 'Client Information', 'Patient Information', and 'VetConnect PLUS'. The 'VetConnect PLUS' tab is active, showing a 'Catalyst Dx Chemistry Analyzer' report. The report includes a table of laboratory results with columns for the test name, value, and reference range. The results are as follows:

Test Name	Value	Reference Range
Creatinine	1.3	0.5 - 1.8 mg/dL
BUN	43	7 - 27 mg/dL
BUN/Creatinine Ratio	33	
Sodium	137	144 - 160 mmol/L
Potassium	10.0	3.5 - 5.8 mmol/L
Chloride	114	109 - 122 mmol/L
ALT	95	10 - 125 U/L

The interface also shows a navigation bar with tabs for 'Test', 'Problems', 'Dr', 'Rx', 'Medical Notes', 'Lab', 'Vaccines', and 'Vital Signs'. A calendar view at the bottom shows the date '5/19/16' and a 'Graphing' option.