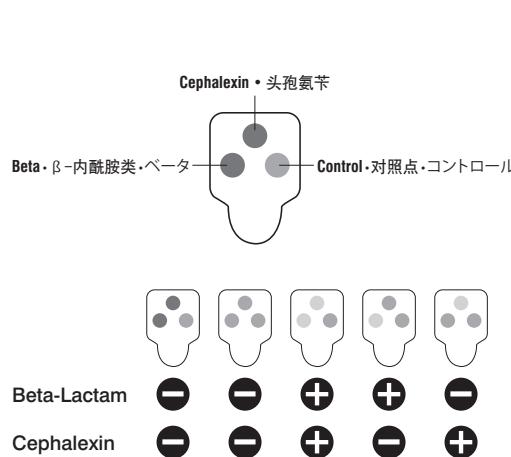
Results

Resultados • Resultados • Résultats • Wyniki • Ergebnisse • Risultati • 结果 • 結果

IDEXX

Simplicity from IDEXX*

It is recommended that positive samples be further analyzed. Cephalexin is a member of the Beta-Lactam class of antibiotics. A positive cephalexin result should be treated as a positive beta-lactam result.

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Technical assistance • idexx.com/dairy

US/Canada/Latinoamerica	+1 800 321 0207	日本	+021 6127 9528
EU	+00800 329 43399	Within AU	+1300 44 33 99
中国	+010 84423133	Within NZ	+0800 102 084

One IDEXX Drive • Westbrook, Maine 04092 USA

*SNAP, SNAPshot and Simplicity from IDEXX are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: idexx.com/patents.

© 2016 IDEXX Laboratories, Inc. All rights reserved. • 06-0004858-01

IDEXX SNAP® Beta-Lactam ST Plus Test

The SNAP® Beta-Lactam ST Plus is a screening test validated to detect beta-lactam residues in commingled cow milk in accordance with European regulations. Review the sensitivity table on the SNAP® Beta-Lactam ST Plus product page on idexx.com/dairy for detection levels of individual drug residues. If you have any questions, please call IDEXX Technical Services. Not for sale in the U.S. Use test before expiration date.

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Storage

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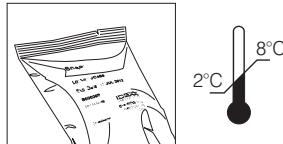
Armazenamento • Stockage

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Ensure test and sample are at correct temperature before use.

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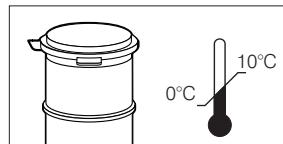
Precautions: Thoroughly mix milk before testing. Do not use abnormal-looking milk or expired tests; do not mix components from different kits. If the control spot fails to develop color, retest the sample.

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1 Transfer milk and swirl until dissolved, but no more than 15 seconds.

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